



## Troubleshooting for Detector baseline


Probable cause	Action
Air bubbles in the inlet side	Tighten connector. Degas solvent.
Worn-out pump seal	Replace pump seal.
Leakage from connector	Retighten leaking part.
Incompletely degassed solvent	Degas solvent again.

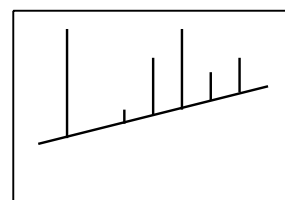


Probable cause	Action
Air bubbles in detector reference	Purge RI detector to replace the reference.
Leakage from connector	Retighten leaking part.

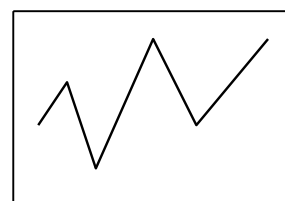

Probable cause	Action
Incompletely degassed solvent	Degas solvent again.

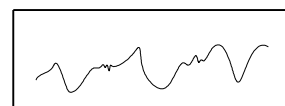

Probable cause	Action
Dust in detector	Clean UV cell. (See section 6.)
For the RI detector, temperature controller failure	Contact our technical service department.

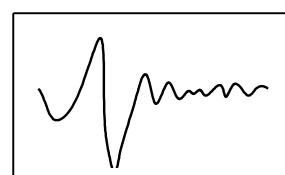

Probable cause	Action
Pump pulsation	Purge the pump.
Unstable solvent flow due to dirt in pump check valve	Disassemble and clean the check valve.

